STOUT™ Recombinant Taq DNA Polymerase Kit

Ordering Info

TBK0022, 20 U (5U/μL)(sample)

TBK0023, 500 U (5U/μL)

TBK0024, 1.000 U (5U/μL)

Description

STOUT™ Recombinant Taq DNA Polymerase Kit is highly flexible kit that includes all PCR components shipped in individual vials to allow for customizable PCR reactions. STOUT Recombinant Taq DNA polymerases (MW 94 kDa) is robust thermostable polymerase produced in *Escherichia coli*. Highly purified the enzyme has a strong 5'→3' polymerase activity. It has a weak 5'→3' exonuclease activity while 3'→5' exonuclease activity is absent.

Features

- **High Thermostability**, the enzyme half-life at 94 °C is 40 minutes.
- **Addition without template** of 3'adenine at the end of PCR fragment.
- **Immediate Activation**.
- Error Rate 1–20x10⁻⁵ errors/bp per cycle.

Applications

- Daily research: standard PCR for targets up 5kb, recombinant clone or cell line checking, etc.
- Generation of PCR fragments for TA cloning.
- DNA Labelling due to capacity to incorporate modified nucleotides.
- DNA Sequencing.

Kit Components

Components	TBK0023	TBK0024
STOUT™ Taq DNA polymerase (5 U/μL)	100 μL	200 μL
STOUT™ PCR Buffer (10x)	1.5 mL	2 x 1.5 mL
MgCl ₂ 25 mM	1.5 mL	2 x 1.5 mL
High-Q™ dNTPs 10 mM TOTAL	1 mL	2 x 1 mL

Order Info Kit Components: STOUT™ Recombinant Taq DNA polymerase (TBZ0201, TBZ0202) | STOUT™ PCR Buffer 10x (TBB0310) | MgCl₂ 25 mM (TBR0215) | High-Q™ dNTPs 10 mM TOTAL (TBR0209).

Storage

Store at -20°C. Shipped in blue ice.

Unit Definition

One unit of STOUT™ Recombinant Taq DNA polymerase catalyzes the incorporation of 10 nanomoles of dNTPs into acid-insoluble material in 30 minutes at 74 °C.

Quality Control

Functionally tested in a 1kb PCR amplification (GC 52%).

Material required (not supplied)

- PCR Grade Water (TBB0303)
- PCR Tubes
- Specific Primers

Also available:

STOUT™ Recombinant Taq DNA Polymerase Master Mix (2x) (TBK0028, TBK0029)

PROTOCOL

1. Thawing all components on ice. Vortex and centrifugate them.
2. On ice, prepare a mix of the following components, considering the number of samples plus two extra reactions.

Reaction Components	Final Concentration	Volume	Volume
STOUT™ PCR Buffer 10x	1 x	2 µL	5 µL
MgCl ₂ 25 mM	2.5 mM	2 µL	5 µL
High-Q™ dNTPs Mix 10 mM _{TOTAL}	0.8 mM	1.6 µL	4 µL
Forward Primer (15 pmol/ µL)	0.2-0.75 µM	0.3-1 µL	1.7-2.5 µL
Reverse Primer (15 pmol/ µL)	0.2-0.75 µM	0.3-1 µL	1.7-2.5 µL
STOUT™ Taq DNA polymerase (5 U/ µL)	0.05 U/ µL	0.2 µL	0.5 µL
Water, molecular biology grade		up 20 µL*	up 50 µL*
DNA template (add in step 4)		*	*
Final Volume		20 µL	50 µL

* consider volume of template to be added in step 4.

3. Distribute the mix prepared in each PCR tube or well.
4. Add in each tube the DNA sample (plasmid DNA: 10-75 ng; gDNA: 100-500 ng). Mix well.
5. Set up thermocycler:

Process	Cycles	Temperature	Time
Initial denaturation	1 x	94 °C	5:00
Denaturation		94 °C	0:35
Annealing	25 - 30 x	T _m	0:35
Extension		72 °C	1:00 per kb
Final Extension	1 x	72 °C	10:00
Conservation	1 x	4 °C	∞

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